


PC BIOTIN ALKYNE

SKU: CCT-1118

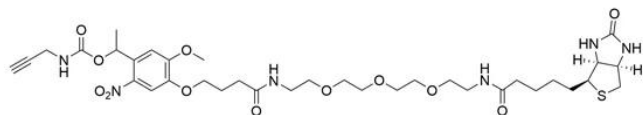

Description

Extraordinary strength of the streptavidin-biotin interaction allows for efficient capturing of even highly dilute targets; however, it makes recovery of proteins from affinity resins challenging. Conventional methods to elute biotinylated proteins from immobilized avidin include the following: (i) denaturation of streptavidin by boiling the resin in a denaturing buffer that may include high concentrations of chaotropic salts, (ii) trypsin digestion of proteins while they are bound to the resin, or (iii) elution of proteins with excess free biotin. These protocols can co-elute contaminant proteins by releasing nonspecifically bound proteins and/or naturally biotinylated proteins concurrently with labeled proteins. In addition, some of these methods can cause elution of high levels of resin-based peptides along with the proteins of interest, resulting in further sample contamination.

PC Biotin Alkyne probes eliminate a major limitation of the streptavidin-biotin affinity purification. This reagent contains a biotin moiety linked to an alkyne moiety through a spacer arm containing a photocleavable linker. Captured biomolecules can be efficiently released under mild, reagent-free conditions (irradiation with near-UV, low intensity lamp) and the small molecular fragment (55.07 Da) left on the labeled protein following cleavage. These features make the photocleavable probe especially attractive for use in biomolecular labeling and proteomic studies.

Specifications

Unit Size	1 mg, 5mg, 25 mg
Molecular weight	780.34
Molecular weight left behind	55.07
Chemical composition	C ₃₅ H ₅₂ N ₆ O ₁₂ S

For research use only. Not intended for animal or human therapeutic or diagnostic use.

CAS	N/A
Solubility	DMSO, DMF, THF, DCM, Chloroform
Appearance	Yellow amorphous solid to yellow oil
Storage Conditions	-20°C.
Shipping Conditions	Ambient temperature

For research use only. Not intended for animal or human therapeutic or diagnostic use.